

# Tuning the Elastic Moduli of Corning® Matrigel® and Collagen I 3D Matrices by Varying the Protein Concentration

CORNING

## Application Note

Katie Slater, Jeff Partridge,  
and Himabindu Nandivada  
Corning Incorporated,  
Life Sciences  
Bedford, MA USA

### Introduction

A complex set of biophysical (e.g., topography, stiffness, viscosity, porosity) and biochemical (e.g., nutrients, matrix composition, cell-matrix interactions) cues control cell biology<sup>1</sup>. In order to mimic the *in vivo* cellular behavior for understanding basic biology, tissue engineering and regenerative medicine applications, it is imperative to better elucidate and control these interactions. The importance of matrix elasticity is being increasingly studied and the stiffness (or elastic modulus) of naturally derived as well as engineered hydrogel substrates has been shown to direct stem cell differentiation<sup>2</sup>, effect cell migration<sup>3</sup>, modulate force of muscle contraction<sup>4</sup>, and influence cell spreading and adhesion<sup>5</sup>. Interestingly, stiffness of tissues in abnormal conditions can also vary: normal rat liver tissue (0.3 to 0.6 kPa) compared to cirrhotic liver (3 to 12 kPa)<sup>6,7</sup>, and normal breast tissue (1.2 kPa) compared to breast tumors (2.4 to 4.8 kPa)<sup>8</sup>. This phenomenon may be causative and not necessarily only an outcome of a disease state<sup>9</sup>.

Lineage specific differentiation of human mesenchymal stem cells (hMSCs) has been shown to vary with matrix stiffness, where hMSCs were directed into particular phenotypes: neurons (matrix stiffness of 0.1 to 1 kPa), myoblasts (8 to 17 kPa), and osteoblasts (25 to 40 kPa)<sup>2</sup>. As rigidity of the cross-linked polyacrylamide gels increased, cells increased their secretory response (e.g., Collagen I). After several weeks of culture and based on matrix rigidity, the specific lineage of the cells was committed. In another study, liver cell differentiation into abnormal phenotypes (myofibroblasts) was driven both by soluble factors and increasing matrix stiffness<sup>6,7</sup>. Following this change, an increased amount of abnormal extracellular matrix (ECM) was produced, resulting in fibrosis (scarring).

Human prostate carcinoma tumor cell (DU-145 cell) migration speed has been influenced by increasing Corning Matrigel matrix stiffness, and the maximum speed was observed at intermediate stiffness<sup>3</sup>. Other factors found to affect migration speed included cell adhesion and proteolytic activity of the cells. Additionally, pore size within the matrix decreased as Matrigel matrix concentration increased; however, this only played a minor role in migration speed.

Increasing Matrigel matrix concentration (from 10% to 40%) was also demonstrated to increase the contraction force of engineered skeletal muscle constructs<sup>4</sup>. Cell survival and function were affected minimally by matrix composition.

### Extracellular Matrices

Basement membrane is a thin layer of ECM proteins underlying all tissues *in vivo*<sup>10</sup>. Corning Matrigel matrix is a solubilized basement membrane preparation extracted from the Engelbreth-Holm-Swarm (EHS) mouse sarcoma, a tumor rich in extracellular matrix proteins, including Laminin (a major component), Collagen IV, heparan sulfate proteoglycans, entactin/nidogen, and a number of growth factors that are found in normal EHS tumors<sup>11,12</sup>. Matrigel matrix has been used in a wide range of applications such as stem cell culture, cell attachment and differentiation, angiogenesis assays, tumor growth, and tissue engineering<sup>13,14</sup>.

Collagen is the most abundant protein present in animals, providing structural and mechanical support<sup>15,16</sup>. There are many different types of Collagen (about 28 known to date)<sup>15</sup> and it is most plentiful in dermis, tendon, and bones<sup>14</sup>. The type I molecule of Collagen is a heterotrimer of 300 nm length and is composed of two  $\alpha_1(I)$  chains and one  $\alpha_2(I)$  chain.

Corning Collagen I rat tail is prepared from rat tail tendons by acid extraction<sup>17</sup>. No enzymes are required for digestion, thus the telopeptides are preserved. Collagen can be used in a thin layer to promote cell attachment, study of tumor cell invasion and migration, fibrillogenesis studies, culture and differentiation of monocytes or macrophages, and in the maintenance of hepatocyte function<sup>18</sup>. Collagen I can also be used in 3D gel form or electrospun into membranes to support 3D cell growth and differentiation<sup>14</sup>.

### **Mechanical Properties of ECMs**

Mechanical properties such as stiffness or elastic modulus of materials can be measured using a variety of techniques such as bulk compressive/tensile testing, rheometry, dynamic light scattering, atomic force microscopy (AFM), and nanoindentation<sup>19,20</sup>.

In rotational or dynamic rheometry, viscoelastic material properties are measured by applying oscillatory shear strain on the sample, most commonly performed in a parallel-plate set-up. For example, in one study, different Corning® Matrigel® matrix (standard product) concentrations ranging from 50% to 100% were tested using a parallel-plate rheometer, and the stiffness ( $G'$ ) linearly increased from approximately 10 to 50 Pa<sup>22</sup>. In a different report, storage modulus measured using a stress-controlled rheometer for Matrigel matrix (high concentration product) concentrations of 4.4, 8, and 17 mg/mL was approximately 20, 70, and 300 Pa, respectively<sup>21</sup>. A stiffness value of 44 Pa has been reported for Matrigel matrix (high concentration product)<sup>22</sup>. Similarly, several publications have reported the stiffness values for Collagen I gels. Using a rotational rheometer, stiffness of Collagen I (rat tail, 2 mg/mL) was reported to be 9 Pa<sup>23</sup>. Furthermore, when Collagen I was mixed with Matrigel matrix at a concentration of either 2 or 4 mg/mL, the stiffness increased to 13.4 and 40.7 Pa, respectively. In another report, the stiffness of Collagen I gels (rat tail and porcine) in the concentration range of 0.5 to 4 mg/mL increased from 0.5 to 100 Pa<sup>24</sup>.

Different mechanical responses are also observed when compressive forces are used instead of shear forces. Using displacement control unconfined compressive testing methodology, a modulus of 6 kPa was achieved for Collagen I (2 mg/mL) samples<sup>25</sup>. The constructs became increasingly stiffer with the addition of Matrigel matrix (10% to 50% addition) with moduli ranging up to 10 kPa. The values obtained in this study are orders of magnitude higher than those from rotational rheometry studies.

Localized mechanical properties can be measured using AFM indentation by pressing a sharp tip into the sample. Under physiological conditions (37°C and aqueous), the average elastic modulus of Matrigel matrix (standard product) measured using AFM with a micron-sized spherical tip was 443 Pa<sup>26</sup>. The differences in the stiffness results between rheological and AFM measurements could be due to differences in length scales (bulk vs. surface) in the two methods.

It should be noted that most of the cell culture work is typically performed on 2D substrates made of rigid materials such as glass and polystyrene, sometimes coated with a thin layer of ECMs. For example, polystyrene and glass substrates have a stiffness value in the GPa range which is orders of magnitude higher than that for ECM proteins (Pa to kPa range) causing the cells to show a non-natural behavior<sup>1</sup>. At the same time, results from measurements of mechanical properties of the thin ECM layers might be confounded by the bulk substrate stiffness.

In this paper, we examined the elastic moduli of the Matrigel matrix and Collagen I gels as a function of the protein concentration. The rheological measurements were performed using single frequency temperature sweeps on a rotational rheometer.

## **Materials and Methods**

### **Corning Matrigel Matrix Preparation**

Corning Matrigel matrix, Growth Factor Reduced (Matrigel GFR, Corning Cat. No. 354230) and Corning Matrigel matrix High Concentration, Growth Factor Reduced (Matrigel HC GFR, Corning Cat. No. 354263) were thawed overnight in ice as directed in the Guidelines for Use<sup>27-29</sup>. The Matrigel matrix products were diluted to the desired protein concentrations using the lot-specific protein concentrations provided on the certificates of analysis using ice-cold Dulbecco's Modified Eagle Medium (DMEM). The contents of the vials were mixed by pipetting up and down and the calculated amount of Matrigel matrix was pipetted using a positive displacement pipet, to ensure accuracy, and added to ice-cold DMEM in pre-chilled glass tubes held in ice. The diluted samples were vortexed gently and held in ice. As polymerization of Matrigel matrix is temperature-dependent and non-reversible, it is important to prepare and hold the samples in ice. The diluted Matrigel matrix samples were mixed by inversion immediately before loading the test samples onto the rheometer using a positive displacement pipet (0.680 mL). Each protein concentration was measured three times.

Samples with concentrations of 3, 6, 9, and 15 mg/mL were made using Matrigel HC GFR. A sample with a concentration of 9 mg/mL was made using Matrigel GFR. The volumes of Matrigel matrix and ice-cold DMEM diluent required were calculated as follows:

$$\frac{\text{Desired final volume (mL)} \times \text{Desired protein concentration (mg/mL)}}{\text{Lot-specific protein concentration (mg/mL)}} = \text{volume of Matrigel matrix required (mL)}$$

$$\text{Desired final volume (mL)} - \text{volume of Matrigel matrix required (mL)} = \text{volume of ice-cold diluent required (mL)}$$

For example, to prepare a 10 mL solution at 9 mg/mL using Corning® Matrigel® matrix HC GFR released with a protein concentration of 19.1 mg/mL:

$$\frac{10 \text{ mL} \times 9 \text{ mg/mL}}{19.1 \text{ mg/mL}} = 4.7 \text{ mL}$$

For example, to prepare a 5 mL solution at 9 mg/mL using Matrigel GFR released with a protein concentration of 9.5 mg/mL:

$$\frac{5 \text{ mL} \times 9 \text{ mg/mL}}{9.5 \text{ mg/mL}} = 4.74 \text{ mL}$$

### Collagen I Preparation

Corning Collagen I High Concentration, rat tail (Collagen HC rat tail, Corning Cat. No. 354249) gelation is pH- and temperature-dependent, and the product must be neutralized in order to polymerize. Collagen HC rat tail was diluted using the alternate gelation procedure included in the lot-specific Certificate of Analysis<sup>30,31</sup>. All dilutions were prepared on ice in pre-chilled glass tubes.

To prepare the desired volume and protein concentration of Collagen I solutions, 0.1X the final volume of 10X Dulbecco's Phosphate Buffered Saline (10X DPBS, Thermo Fisher), 0.023 mL of 1N Sodium Hydroxide (JT Baker) per mL of Collagen I were added and the final volume was adjusted using ice-cold DI-water. The contents of the tube were mixed by vortexing and the required volume of Collagen I was added to the solution using a positive displacement pipet, mixed by vortexing gently and held on ice. All diluted samples were mixed by inversion immediately before loading the test samples onto the rheometer using a positive displacement pipet (0.680 mL). Each protein concentration was measured in triplicate.

Samples with concentrations of 1, 3, and 7 mg/mL were made using Collagen HC rat tail.

For example, to prepare a 6 mL solution with a final Collagen I concentration of 7 mg/mL using Collagen HC rat tail (Lot No. 7016285) with a protein concentration of 9.82 mg/mL:

*0.6 mL of 10X DPBS, 0.0984 mL of 1N NaOH and 1.02 mL of ice-cold DI-water were added to a pre-chilled glass tube and mixed by vortexing. Using a positive displacement pipet, 4.28 mL of Collagen HC rat tail was added to the glass tube.*

### Rheological Measurements

The samples were analyzed using the Kinexus® Ultra+ Rotational Rheometer (Malvern Instruments) using a parallel plate (40 mm roughened) configuration. Using a positive displacement pipet, samples (0.680 mL) were directly added to center of the lower plate at a temperature of 5°C. The top plate was immediately lowered before the gel started to form, to a working gap of 0.5 mm. The temperature was ramped from 5°C to 37°C at a rate of 10°C/minute and the temperature was held at 37°C. A solvent trap was used to cover the sample and prevent any evaporation effects around the perimeter. Viscoelastic measurements were performed periodically at a single frequency of 0.5 Hz and 0.2% strain. Measurements were recorded until the elastic modulus reached equilibrium. The data was acquired and analyzed using the rSpace® software (Malvern Instruments).

### *Tips for Successful Usage of Corning® Matrigel® Matrix*

#### **Storage**

- ▶ Store in a -20°C non-frost-free freezer.
- ▶ Minimize temperature fluctuations, do not store on the door or in a freezer that is opened frequently.

#### **Thawing**

- ▶ Pack in ice (completely submerge).
- ▶ Cover ice bucket and place in 2°C to 8°C location (cold room or refrigerator) overnight.
- ▶ Avoid repeated freeze thaw cycles.
- ▶ Thaw one-time use aliquots using the same guidelines, time may be adjusted based on the volume of the aliquot.

#### **Diluting**

- ▶ Use lot-specific protein concentration provided on the Certificate of Analysis to prepare dilutions.
- ▶ Keep Matrigel matrix and all solutions on ice.
- ▶ Add Matrigel matrix to ice-cold solutions (do not dilute with water, aggregates may form).
- ▶ Use pre-chilled tips, tubes, and cultureware.
- ▶ Use positive displacement pipets for accurate pipetting (due to viscosity).

#### **Handling**

- ▶ Plan ahead.
- ▶ Thaw Matrigel matrix appropriately.
- ▶ Make one-time use aliquots.
- ▶ Prepare Matrigel matrix aliquots immediately after thawing the vial.
- ▶ Store one-time use aliquots using the same recommended storage conditions.
- ▶ Aliquot non-diluted Matrigel matrix only.
- ▶ Matrigel matrix is temperature-sensitive and will start to gel above 10°C.
- ▶ Polymerization is non-reversible.

### *Tips for Successful Usage of Corning Collagen I, Rat Tail*

#### **Storage**

- ▶ Store at 2°C to 8°C.
- ▶ Do not freeze the product.

#### **Diluting**

- ▶ Use lot-specific protein concentration provided on the Certificate of Analysis to prepare dilutions.
- ▶ Use ice-cold solutions.
- ▶ Keep product and diluted samples on ice.
- ▶ Use positive displacement pipets for accurate pipetting (due to viscosity).


#### **Handling**

- ▶ Polymerization is pH- and temperature-dependent.
- ▶ Collagen must be neutralized in order to form a gel.

## **Results and Discussion**


Proteins present in naturally derived ECMs such as Corning Matrigel matrix and Collagen I self-assemble into complex 3D fibrous structures<sup>32</sup>. Mechanical properties of these matrices regulate a myriad of cellular functions including cell spreading, proliferation, migration, and differentiation<sup>11</sup>. Mechanical properties such as stiffness or elastic modulus of materials can be measured using a variety of techniques such as bulk compressive/tensile testing, rheometry, AFM, and nanoindentation.

To investigate whether varying the protein concentration alters the elastic modulus of the Matrigel matrix gels, Corning Matrigel High Concentration, Growth Factor Reduced (Matrigel HC GFR; at a concentration of 19.1 mg/mL) was diluted using ice-cold DMEM and liquid samples were directly added to the rotational rheometer. The temperature of the bottom plate (Peltier-controlled) was ramped from 5°C to 37°C and then held at 37°C, which is the recommended gelation temperature for Matrigel matrix. The evolution of the elastic component ( $G'$ ) and the viscous component ( $G''$ ) of the complex shear modulus were observed over time and the sample was allowed to reach a state of equilibrium. Over a range of concentrations from 3 to 19.1 mg/mL for Matrigel HC GFR, the elastic component of the shear modulus increased from  $9.1 \pm 0.3$  Pa to  $288.2 \pm 9$  Pa (Figures 1A and 1B). At the equilibrium state, the plateau value of the elastic component was larger than the viscous component indicating that the gel finally formed was elastic in nature (Figure 1C).

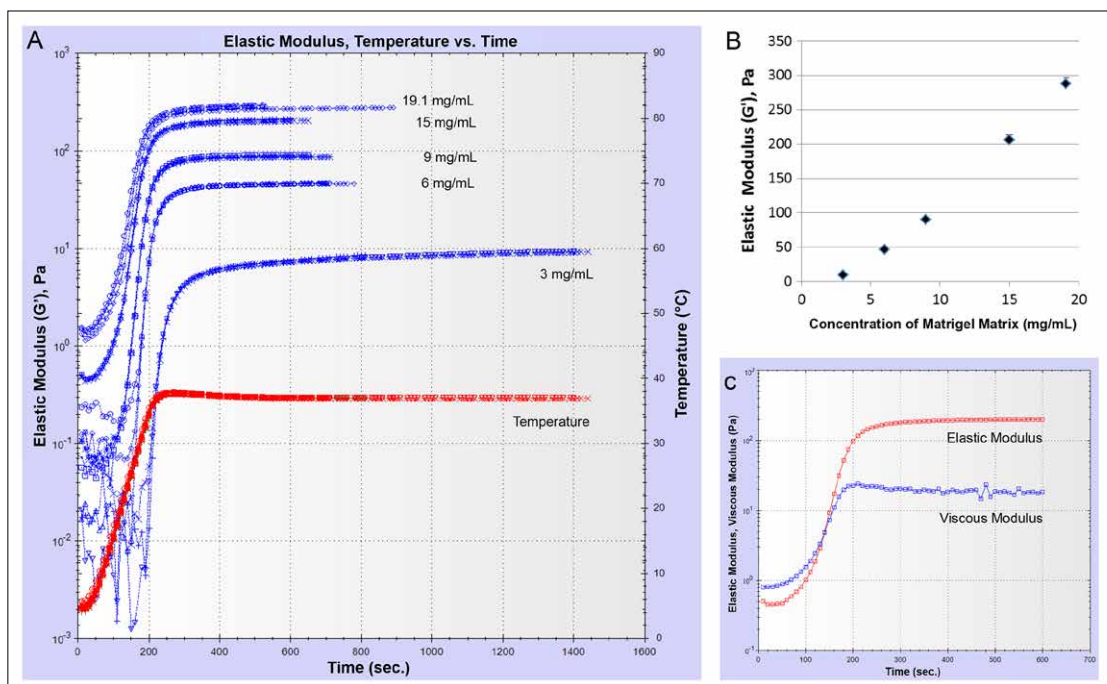

**Figure 1. Increase of elastic modulus as a function of Corning Matrigel HC GFR concentration.** (A) A plot of the elastic modulus ( $G'$ ) as a function of time for samples of Matrigel matrix at different concentrations. Temperature was ramped from 5 $^{\circ}\text{C}$  to 37 $^{\circ}\text{C}$  and held at 37 $^{\circ}\text{C}$ . (B) Plateau values of  $G'$  (in Pa) shown as a function of Matrigel matrix concentration ( $n = 3$ ). (C) Representative time course of elastic modulus and viscous modulus for Matrigel matrix at a concentration of 15 mg/mL.


Corning<sup>®</sup> Matrigel<sup>®</sup> matrix is available as Matrigel HC GFR (a high concentration product) with a concentration range of 18 to 22 mg/mL and Corning Matrigel Growth Factor Reduced (Matrigel GFR) with a concentration range of 7 to 10 mg/mL. The shear moduli (elastic component) of gels formed using the Matrigel HC GFR and Matrigel GFR at the same protein concentration (9 mg/mL) were measured and found to be comparable (Figure 2) indicating that mechanical properties of Matrigel matrix are not influenced by the small differences in the processing of the high concentration and standard concentration products.

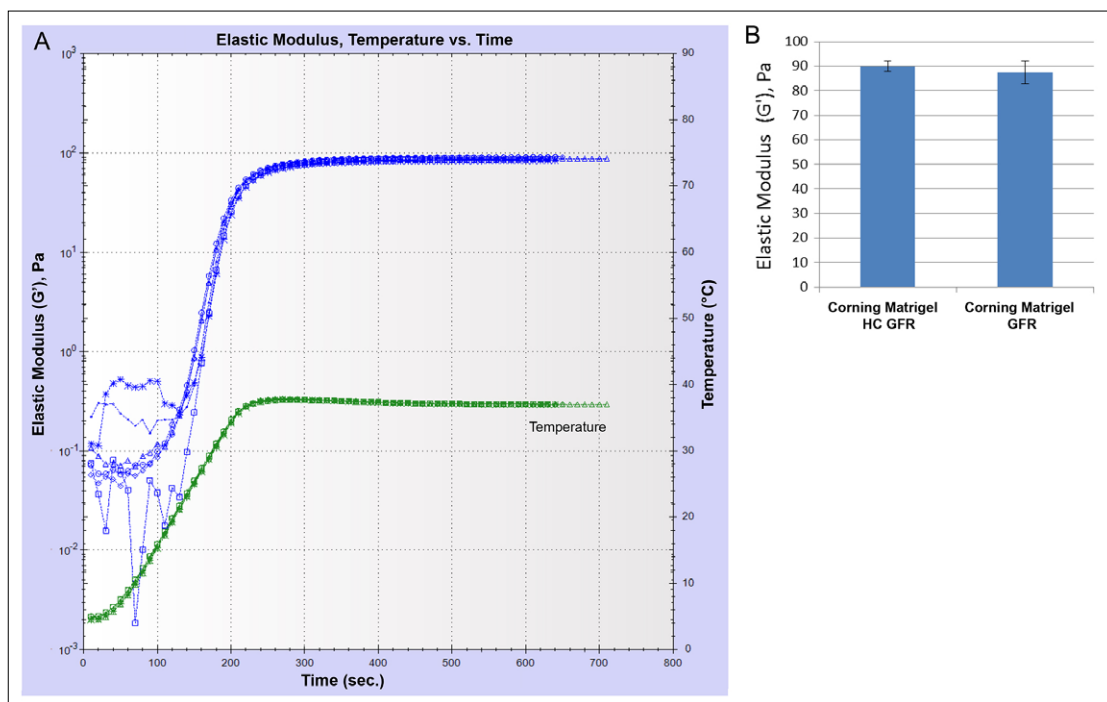

**Figure 2. Comparison of elastic modulus ( $G'$ ) for two different Corning Matrigel matrix products formed at the same protein concentration (9 mg/mL).** (A) A plot of the elastic modulus as a function of time for samples of Matrigel HC GFR and Matrigel GFR. Temperature was ramped from 5 $^{\circ}\text{C}$  to 37 $^{\circ}\text{C}$  and held at 37 $^{\circ}\text{C}$ . (B) Plateau values of  $G'$  (in Pa) shown for the two different Matrigel matrix products ( $n = 3$ ).

To investigate whether varying the protein concentration alters the elastic modulus of the Collagen I gels, samples of different concentrations were created by neutralizing Corning® Collagen I High Concentration, rat tail (Collagen HC rat tail; at a concentration of 9.82 mg/mL) with 1 N NaOH solution. For Collagen HC rat tail concentrations of 1, 3, and 7 mg/mL,  $G'$  plateaued at  $5.0 \pm 0.6$  Pa,  $55.4 \pm 11.0$  Pa, and  $341.8 \pm 32.4$  Pa, respectively (Figures 3A and 3B). Collagen I gels demonstrated an elastic behavior similar to Matrigel matrix gels as seen when the plateau values of  $G'$  and  $G''$  were compared (Figure 3C).

Other factors that also play an important role in the gel formation and thereby affect the modulus are gelation temperature, temperature ramp, changes in general handling of the protein solutions, and more specifically for Collagen I pH of the solution. Rheometer parameters such as applied shear, frequency, working gap, and also the type of instrument used for the measurement could affect the final modulus values attained.

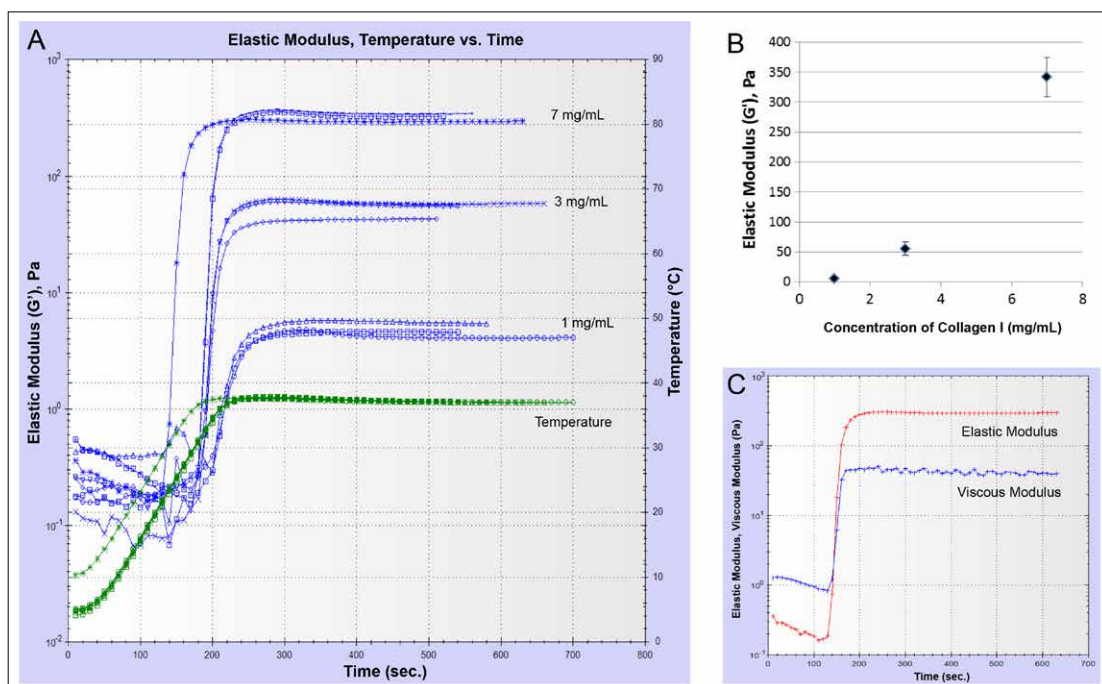

**Figure 3. Increase of elastic modulus as a function of Collagen HC rat tail concentration.** (A) A plot of the elastic modulus ( $G'$ ) over time for samples of Collagen I at different concentrations. Temperature was ramped from 5°C to 37°C and held at 37°C. (B) Plateau values of the shear modulus (elastic component,  $G'$ ) (Pa) vs. Collagen I concentration ( $n = 3$ ). (C) Representative time course of elastic modulus and viscous modulus for Collagen I at a concentration of 7 mg/mL.

## Conclusions

Cellular behavior is governed by a complex interplay of various physico-chemical factors. For instance, the effects of mechanical properties of substrates on cellular function with respect to understanding cell biology, and in applications of cell therapy and regenerative medicine continue to be an area of intense study. Numerous highly sensitive tools and techniques are available for the measurement of the mechanical properties of matrices. In this paper, we investigated the effect of protein concentration on the elastic moduli of Corning Matrigel® matrix and Collagen I gels using a rotational rheometer. It was observed that the elastic modulus of both materials increased with increasing protein concentration and the gels were elastic in nature.

## Acknowledgments

We are grateful to Dr. Philip Rolfe and Maria Lightfoot (Malvern Instruments, Westborough, MA) for giving us access to the Kinexus® Ultra+ Rheometer and assisting with set-up, data collection, and analysis.

## References

1. Santos, E. et al. Cell-biomaterial interaction: Strategies to mimic the extracellular matrix. Chapter 25, on biomimetics, Pramatarova L. (Ed.), InTech (2011).
2. Engler, A. J. et al. Matrix elasticity directs stem cell lineage specification. *Cell*, 126, 677-689 (2006).
3. Zaman, M. H. et al. Migration of tumor cells in 3D matrices is governed by matrix stiffness along with cell-matrix adhesion and proteolysis. *Proceedings of National Academy of Sciences USA*, 103(37) (2006).
4. Hinds, S. et al. The role of extracellular matrix composition in structure and function of bioengineered skeletal muscle. *Biomaterials*, 32, 3575-3583 (2011).
5. Pelham, Jr R. J. and Wang Y-L. Cell locomotion and focal adhesions are regulated by substrate flexibility, *Proceedings of National Academy of Sciences USA*, 94, 13661-13665, (1997).
6. Olsen, A. L. et al. Hepatic stellate cells require a stiff environment for myofibroblastic differentiation, *American Journal of Physiology - Gastrointestinal and Liver Physiology*, 301, G110-G118 (2011).
7. Li, Z. et al. Transforming growth factor- $\beta$  and substrate stiffness regulate portal fibroblast activation in culture, wells. *Hepatology*, 46(4), 1246-1256 (2007).
8. Keely, P. and Nain, A. Capturing relevant extracellular matrices for investigating cell migration. *F1000Research*, 4 (F1000 Faculty Rev), 1408 (2005).
9. Liu, F. et al. Feedback amplification of fibrosis through matrix stiffening and COX-2 suppression. *Journal of Cell Biology*, 190(4), 693-706 (2010).
10. LeBleu, V. S. et al. Structure and function of basement membranes. *Experimental Biology and Medicine*, 232(9), 1121-1129 (2009).
11. Kleinman, H. K. et al. Isolation and characterization of type IV procollagen, laminin, and heparan sulfate proteoglycan from the EHS sarcoma. *Biochemistry*, 21, 6188 (1982).
12. Kleinman, H. K. et al. Basement membrane complexes with biological activity. *Biochemistry*, 25, 312 (1986).
13. Kleinman, H. K. and Martin, G. R. Matrigel: Basement membrane matrix with biological activity. *Seminars in Cancer Biology*, 378-386 (2005).
14. Sanyal, S. Culture and assay systems used for 3D cell culture. Corning CLS-AC-AN-245.
15. Shoulders, M. D. and Raines, R. T. Collagen structure and stability. *Annual Review of Biochemistry*, 78, 929-958 (2009).
16. Muiznieks, L. D. and Keeley, F. W. Molecular assembly and mechanical properties of the extracellular matrix: A fibrous protein perspective. *Biochimica et Biophysica Acta- Molecular Basis of Disease*, 1832, 866-875 (2013).
17. Artym, V. V. and Matsumoto, K. Imaging cells in three-dimensional collagen matrix. *Current Protocols in Cell Biology*, 48:10.18, 10.18.1-10.18.20 (2010).
18. Cheema, U. et al. Collagen: Applications of a natural polymer in regenerative medicine. *Regenerative Medicine and Tissue Engineering - Cells and Biomaterials*, Eberli D. (Ed.), InTech (2011).
19. Vinckier, A. and Semenza, G. Measuring elasticity of biological materials by atomic force microscopy. *FEBS Letters*, 430(1-2), 12-16 (1998).
20. McKee, C. T. et al. Indentation versus tensile measurements of Young's modulus for soft biological tissues. *Tissue Engineering Part B*, 17(3), 155-164 (2011).
21. Chaudhuri, O. et al. Extracellular matrix stiffness and composition jointly regulate the induction of malignant phenotypes in mammary epithelium. *Nature Materials*, 13, 970-978 (2014).
22. Lambrecht, L. et al. The type and composition of alginate and hyaluronic-based hydrogels influence the viability of stem cells of the apical papilla. *Dental Materials*, 30(12):e349-61 (2014).
23. Anguiano, M. et al. Characterization of three-dimensional cancer cell migration in mixed collagen-Matrigel scaffolds using microfluidics and image analysis. *PLoS ONE*, 12(2): e0171417 (2017).
24. Motte, S. and Kaufman, L. J. Strain stiffening in Collagen-I networks, *Biopolymers*, 99(1), 35-46 (2012).
25. Dewitt, D. D. et al. Collagen I-Matrigel scaffolds for enhanced schwann cell survival and control of three-dimensional cell morphology. *Tissue Engineering Part A*, 15(10), 2785-2793 (2009).
26. Soofi, S. S., et al. The elastic modulus of Matrigel as determined by atomic force microscopy. *Journal of Structural Biology*, 167(3), 216-219 (2009).
27. Corning Matrigel Basement Membrane Matrix Growth Factor Reduced, 10 mL; Guidelines for Use (SPC-354230).
28. Corning Matrigel Basement Membrane Matrix High Concentration Growth Factor Reduced, 10 mL; Guidelines for Use (SPC-354263).
29. Corning Matrigel Matrix Frequently Asked Questions (CLS-DL-CC-026).
30. Corning Collagen I High Concentration (HC), Rat Tail; Certificate of Analysis (SPC-354249).
31. Corning Collagen Products Frequently Asked Questions (CLS-DL-AC-004).
32. Trappmann, B. and Chen, C. How cells sense extracellular matrix stiffness: a material's perspective. *Current Opinion in Biotechnology*, 24(5), 948-953 (2013).

For more specific information on claims, visit the Certificates page at [www.corning.com/lifesciences](http://www.corning.com/lifesciences).

**Warranty/Disclaimer:** Unless otherwise specified, all products are for research use only. Not intended for use in diagnostic or therapeutic procedures. Not for use in humans. Corning Life Sciences makes no claims regarding the performance of these products for clinical or diagnostic applications.

For additional product or technical information, visit [www.corning.com/lifesciences](http://www.corning.com/lifesciences) or call 800.492.1110. Outside the United States, call +1.978.442.2200 or contact your local Corning sales office.

# CORNING

## Corning Incorporated Life Sciences

836 North St.  
Building 300, Suite 3401  
Tewksbury, MA 01876  
t 800.492.1110  
t 978.442.2200  
f 978.442.2476

[www.corning.com/lifesciences](http://www.corning.com/lifesciences)

### ASIA/PACIFIC

**Australia/New Zealand**  
t 61 427286832

**China**  
t 86 21 3338 4338  
f 86 21 3338 4300

**India**  
t 91 124 4604000  
f 91 124 4604099

### Japan

t 81 3-3586 1996  
f 81 3-3586 1291

### Korea

t 82 2-796-9500  
f 82 2-796-9300

### Singapore

t 65 6572-9740  
f 65 6735-2913

### Taiwan

t 886 2-2716-0338  
f 886 2-2516-7500

### EUROPE

CSEurope@corning.com

### France

t 0800 916 882  
f 0800 918 636

### Germany

t 0800 101 1153  
f 0800 101 2427

### The Netherlands

t 020 655 79 28  
f 020 659 76 73

### United Kingdom

t 0800 376 8660  
f 0800 279 1117

### All Other European Countries

t +31 (0) 206 59 60 51  
f +31 (0) 206 59 76 73

### LATIN AMERICA

grupoLA@corning.com

### Brasil

t 55 (11) 3089-7400

### Mexico

t (52-81) 8158-8400