

CORNING

# Pipet Controller Charging Stand

For Corning® Stripettor® Ultra and Pro Pipet Controllers


**Instruction Manual**

Catalog Number:  
4998




## 1. Introduction

The Pipet Controller Charging Stand for Corning® Stripettor® Ultra and Pro Pipet Controllers is intended to conveniently charge and store a pipet controller when not in use. Depending on the use needs, there are three positions to store the pipet controller, and are visible as icons printed on the stand.

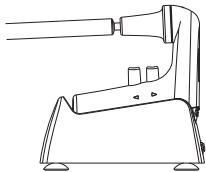


### **Break position**

Convenient while working, helps to avoid contamination.

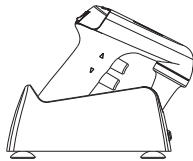

### **Rest position**

Ideal way to store a pipet controller after use.

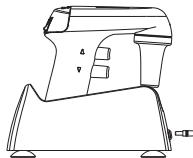

### **Charging position**

Used while pipet controller is in charging process.

## 2. Packaging

- ▶ 3-position charging stand
- ▶ Instruction manual


## 3. Charger

### IMPORTANT:

- ▶ The pipet controller may be charged only with the original charger.
- ▶ The mains voltage shall conform with the charger specification (Input: 100-240V, 50/60Hz, 0.2A; Output: 9V).

## 4. Charging the batteries on the charging stand

- ▶ Put the pipet controller on the stand.
- ▶ Plug the charger into the stand outlet.
- ▶ Connect the charger to the adapter and to the main power.

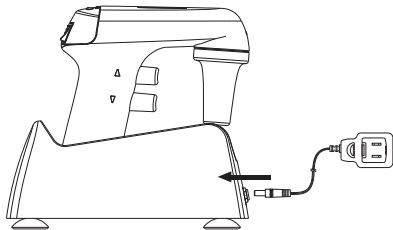

## 5. Equipment Disposal

According to Directive 2012/19/EU of the European Parliament and Council of 4th July 2012 on waste electrical and electronic equipment (WEEE) as amended, this product is marked with the crossed-out wheeled bin and must not be disposed of with domestic waste.

Consequently, the buyer shall follow the instructions for reuse and recycling of waste electronic and electrical equipment (WEEE) provided with the products and available at the following link: **[www.corning.com/weee](http://www.corning.com/weee)**.

For more information about the charging process visit **[www.corning.com/lifesciences](http://www.corning.com/lifesciences)** or user manuals of respective products.

**CORNING**

**Corning Incorporated**

*Life Sciences*

**[www.corning.com/lifesciences](http://www.corning.com/lifesciences)**