

DMSO Removal and CD3⁺ Cell Selection from Cryopreserved Apheresis **Products Using Novel BACS Technology**

Content originally prepared for ThermoGenesis by Dalip Sethi¹, Jillian Miller¹, Zelenia Contreras², John Perea², and Stephen Truong² ¹Cesca Therapeutics, Inc., Rancho Cordova, CA USA; ²ThermoGenesis Corp., Rancho Cordova, CA USA

Introduction


Cancer therapy is a complex and growing field of medicine. With the advent of adoptive immunotherapy, coupled with excellent clinical outcomes in CD19⁺ hematologic malignancies, CAR-T cell therapy has created intense interest within the medical and scientific communities. Although genetic-modification of immune cells is one of the most advanced therapies, the manufacturing process still relies on traditional cell processing techniques. We have previously developed a guick and easy method to remove DMSO from thawed apheresis products and have now coupled this technology with the selection of CD3⁺ T-cells using ThermoGenesis' technology, Buoyancy-Activated Cell Sorting (X-BACS[™]) process. DMSO removal was successfully achieved using the X-WASH® System.

Materials and Methods


Healthy adult cryopreserved apheresis samples were thawed at 37°C and diluted with thaw-wash media (2.5% HSA, 5% Dextran, and DNase I in saline). The sample was transferred to an X-WASH Disposable Cartridge. Post-centrifugation, waste media was removed using positive pressure, and new media was added using controlled negative pressure. Washed cells were transferred to a new Disposable Cartridge and incubated with anti-CD3 antibody followed by incubation with microbubble reagent. The target and non-target fractions were separated using centrifugation. CD3+ cells were also studied for expansion using CD3/CD28 and IL-2 activation.

In a separate experiment, the expansion of CD3⁺ cells isolated using X-BACS was equivalent to CD3⁺ cells isolated using technologies from other manufacturers. Viable CD3⁺ cells were quantified over 14 days using flow cytometry.

CORNING

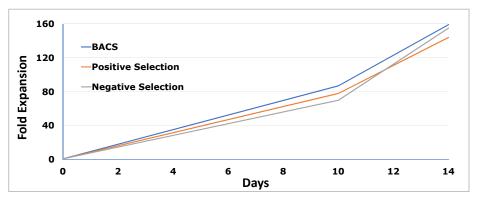


To assess the consistency of CD3⁺ cell recoveries using the X-BACS process, the antibody and microbubble concentrations were kept constant while cell concentrations were increased. The results indicate only a marginal drop in CD3⁺ cell recovery within the normally expected dose range (30 to 200 x 10⁶/mL).

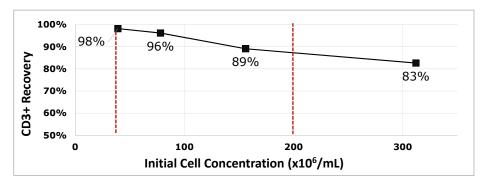

Conclusions


DMSO removal using the X-WASH System resulted in cell recoveries of greater than 90% with no significant loss of viability. A mean CD3⁺ cell recovery of 84% was accomplished using the X-BACS process with high purity (>94%). In brief, we have developed an efficient method for removing DMSO and isolating CD3⁺ T-cells from cryopreserved apheresis samples that can be used for CAR-T cell manufacturing.

Process Flow

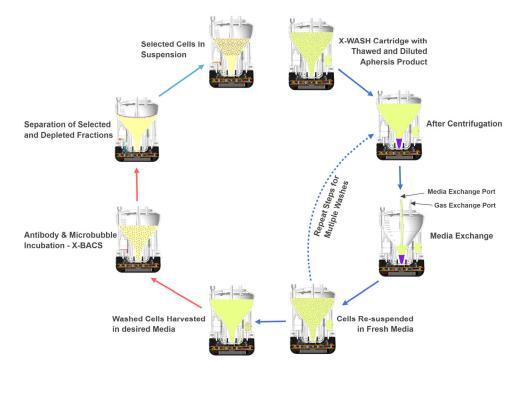

Results and Discussion

The DMSO removal from cryopreserved samples and CD3⁺ cell isolation was accomplished in two phases. Each phase was studied for recovery efficiency and cell viability.

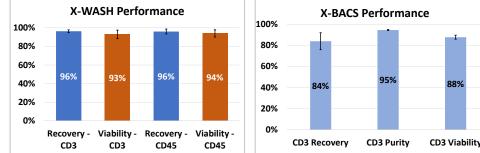

Overall recovery was calculated by multiplying values obtained in X-WASH and X-BACS steps. A mean CD3⁺ cell recovery of 81% was observed (n=3).

	Overall Performance Data	
CD3 Viability	88%	
CD3 Purity	95%	I
CD3 Recovery	81%	

that made by Corning the void and a control and medical over a support that claim and that you want to be the product of any medicad action and that you want to be the product of any medicad action and the product of any medicad action and that you want to be the product of any medicad action and the product of any medicad action action