Corning[®] rLaminin-521 (Human)

CORNING

Corning has partnered with BioLamina of Sweden for the supply of recombinant human laminin-521. Corning rLaminin-521 (Human) is a heterotrimer composed of α 5, β 2, and γ 1 chains expressed in a mammalian cell culture system. rLaminin-521 (Human) supports long-term self-renewal of human pluripotent stem cells (hPSC), including embryonic stem cells (hESC) and induced pluripotent stem cells (iPSC) in defined and xeno-free environments. rLaminin-521 provides additional benefits, including ROCK inhibitor independent single cell expansion of PSCs, which improves hPSC culture ease and efficiency.

Features

- Recombinant protein
- Easy to self-coat
- Available in 0.1 mg aliquots for easy handling
- > Compatible with multiple hPSC culture media and dissociation reagents
- Supports undifferentiated expansion of hPSC
- > Suitable for xeno-free, defined culture environment



Human Pluripotent Stem Cell Culture (hPSC)

hPSC, including embryonic stem cells and induced pluripotent cells, exhibit unlimited self-renewal capability and differentiation potential. These cells can give rise to the majority of cell types in the body and can be a source of cells and tissues for regenerative medicine. hPSC-derived cells can also be used as a model for drug screening. Typically feeder-free hPSC culture is performed using clump passage, as single cell culture requires the addition of ROCK inhibitor in the media. Long term effects of using these chemicals are unknown. Clump passage methodology is uncontrolled and not convenient to scale-up.

Morphology

hESC cultured on rLaminin-521 (Human) in xeno-free medium exhibited characteristic colony morphology with a high nuclear-to-cytoplasm ratio.





20x magnification

10x magnification

Marker Expression

Markers for undifferentiated hPSC were determined by immunocytochemistry. Image analysis indicates Oct-4 (green) expression in the cells. Nuclei were stained with DAPI (blue).



Oct-4 + DAPI

Tri-Lineage Differentiation

After 10 passages in xeno-free medium, pluripotency of hESC were assessed by differentiation into embryoid bodies (EBs) and quantification of lineage-specific markers in EBs by quantitative RT PCR. Markers from all 3 lineages were expressed in EBs, which demonstrated that the cells maintained pluripotency after multiple passages.



Karyotype Analysis

hESC cultured on rLaminin-521 (Human) in xeno-free medium maintained normal karyotype. Karyotyping was performed at multiple times during the expansion of the cells.



hPSC Lines, Media, and Dissociation Reagents Evaluated on rLaminin-521 (Human)

Cells	Human Embryonic Stem Cells, Human Induced Pluripotent Stem Cells		
Media	mTeSR™1², TeSR™2², Essential E8¹, NutriStem™ XF/FF¹,2		
Dissociation reagents	TrypLE [™] Select ² , EDTA ¹ , Trypsin/EDTA ²		

¹ Lu, et al., (2014). A defined xeno-free and feeder-free culture system for the derivation, expansion and direct differentiation of transgene-free patient-specific induced pluripotent stem cells. Biomaterials 35:2816.

² Rodin, et al., (2014). Clonal culturing of human embryonic stem cells on laminin-521/E-cadherin matrix in defined and xeno-free environment. Nature Communications 5:3195.

Ordering Information

Corning® rLaminin-521 (Human)

Cat. No.	Description	Qty/Pk	Qty/Cs
354220	20 μg vial	1	1
354221	100 μg vial	1	1
354222	1 mg (10 x 100 μg vials)	1	1
354223	5 mg (50 x 100 μg vials)	1	1

Warranty/Disclaimer: Unless otherwise specified, all products are for research use only. Not intended for use in diagnostic or therapeutic procedures. Not for use in humans. Corning Life Sciences makes no claims regarding the performance of these products for clinical or diagnostic applications.



At Corning, cells are in our culture. In our continuous efforts to improve efficiencies and develop new tools and technologies for life science researchers, we have scientists working in Corning R&D labs across the globe, doing what you do every day. From seeding starter cultures to expanding cells for assays, our technical experts understand your challenges and your increased need for more reliable cells and cellular material.

It is this expertise, plus a 160-year history of Corning innovation and manufacturing excellence, that puts us in a unique position to offer a beginning-to-end portfolio of high-quality, reliable cell culture consumables.

www.corning.com/lifesciences/solutions

For additional product or technical information, please visit www.corning.com/lifesciences or call 800.492.1110. Customers outside the United States, please call +1.978.442.2200 or contact your local Corning sales office.

Corning Incorporated Life Sciences 836 North St. Building 300. Suite 3401	Worldwide Support Offices ASIA/PACIFIC	Japan t 81 3-3586 1996 f 81 3-3586 1291 Korea	EUROPE France t 0800 916 882 f 0800 918 636	All Other European Countries t 31 (0) 20 659 60 51 f 31 (0) 20 659 76 73
Tewksbury, MA 01876 t 800.492.1110	Australia/New Zealand t 0402-794-347	t 82 2-796-9500 f 82 2-796-9300	Germany t 0800 101 1153 f 0800 101 2427 The Netherlands t 31 20 655 79 28	LATIN AMERICA Brasil t (55-11) 3089-7419 f (55-11) 3167-0700 Mexico t (52-81) 8158-8400 f (52-81) 8313-8589
t 978.442.2200	China t 86 21 2215 2888 f 86 21 6215 2988	Singapore t 65 6733-6511 f 65 6861-2913		
f 978.442.2476 www.corning.com/lifesciences				
	India t 91 124 4604000 f 91 124 4604099	Taiwan t 886 2-2716-0338 f 886 2-2516-7500	f 31 20 659 76 73	
			United Kingdom t 0800 376 8660 f 0800 279 1117	

GOSSELIN

PYREX

FALCON

Laminin-521 is manufactured by BioLamina AB, Sweden for Corning. For a listing of trademarks, visit us at www.corning.com/lifesciences/trademarks.

CORNING |

All other trademarks are the property of their respective owners.