

# Corning® Fresh Human Hepatocyte Cytochrome P450 Induction Assays

CORNING

## Protocol - Hepatocyte Assay Methods

The cellular content of cytochromes P450 can be induced several fold by pre-treatment with drugs and other xenobiotics. Drug-mediated induction of P450 can result in increased metabolism of other drugs or itself, leading to potentially harmful drug interactions and/or drug tolerance. Freshly isolated human hepatocytes provide an excellent *in vitro* means for predicting P450 induction potential by drug candidates.

### Purpose

The following protocol describes a method for treating plated human hepatocytes with “classical” P450 inducers (i.e. Rifampicin-CYP3A4, 2C9/19; Phenobarbital-CYP2B6, 3A4; and  $\beta$ -Naphthoflavone-CYP1A2) and analysis of results. The same general protocol can be followed when testing the induction potential of new chemical entities (NCEs).

**Table 1: Chemicals and Suppliers**

Chemical Reagent	Supplier (Cat. No.)
Rifampicin (RIF)	Sigma Chemical Co. (R-3501)
$\beta$ -naphthoflavone ( $\beta$ -NF)	Sigma Chemical Co. (N-3633)
Phenobarbital (PB)	Sigma Chemical Co. (P-5178)
Corning Hepatocyte Culture Media Kit	Corning Life Sciences (355056)
Dimethyl Sulfoxide (DMSO)	JT Baker (T15599)
KHB Buffer	Sigma Chemical Co. (K-3753)
Phenacetin (CYP1A2 substrate)	Sigma Chemical Co. (A-2375)
Testosterone (CYP3A4 substrate)	Sigma Chemical Co. (T-1500)
Diclofenac (CYP2C9 substrate)	Sigma Chemical Co. (D-6899)
[14C]S-Mephenytoin (CYP2C19/2B6 substrate)	GE Life Sciences (CFA-763)
S-Mephenytoin (CYP2C19/2B6 substrate)	Corning Life Sciences (451032)

**Table 2: P450 Inducers and Substrates/Solvents and Concentrations**

Chemical	Stock Concentration	Solvent	Final Working Concentration
Rifampicin (RIF)	25 mM	DMSO	20 $\mu$ M
$\beta$ -Naphthoflavone ( $\beta$ -NF)	25 mM	DMSO	20 $\mu$ M
Phenobarbital (PB)	0.2 M	PBS	2 mM
Testosterone	0.1 M	DMSO	200 $\mu$ M
Diclofenac	0.1 M	DMSO	100 $\mu$ M
Phenacetin	0.1 M	DMSO	100 $\mu$ M
*[14C]S-Mephenytoin	10 mM	ACN	100 $\mu$ M

\* Specific Activity of 10 mM stock: 5 to 7 mCi/mmol.

### Other Material and Equipment:

- ▶ Biosafety hood
- ▶ Incubator with 37°C and 95% air/5% CO<sub>2</sub> at atmospheric level capacity
- ▶ Basic cell culture equipment

### Fresh Plated Hepatocytes:

#### Plated Human Hepatocytes/Corning® BioCoat™ Collagen I microplates

- ▶ 6 well Corning BioCoat Collagen I plates (Cat. No. 454406)
- ▶ 12 well Corning BioCoat Collagen I plates (Cat. No. 454412)
- ▶ 24 well Corning BioCoat Collagen I plates (Cat. No. 454424)
- ▶ 48 well Corning BioCoat Collagen I plates (Cat. No. 454425)
- ▶ 96 well Corning BioCoat Collagen I plates (Cat. No. 454496)

### Procedure

1. Unpack the plated hepatocytes. Follow enclosed instructions for handling requirements and maintenance until ready for further testing.
2. Incubate the plated hepatocytes for 1-2 days in a humidified 37°C incubator with 95% air/5% CO<sub>2</sub> atmosphere to allow for recovery from shipping.
3. Prepare the chemical inducers (RIF, β-NF, and PB), solvent vehicle controls (DMSO for RIF and β-NF, and PBS for PB) and any test compounds (Table 3). The media/inducer preparation should be prepared fresh on the day of use.

**Note:** Whenever DMSO is used to dissolve test compound the final DMSO concentration should not exceed 0.1%.

**Table 3: Treatment Schedule and Inducer Concentrations**

Chemical Inducer	Working Concentration	P450	Total Treatment Time
Rifampicin	20 μM	3A4, 2C9, 2C19	72 hours
PB	2 mM	2B6	72 hours
β-NF	20 μM	1A2	72 hours

4. Aspirate the media from the well and replace the media containing the desired treatment condition. Use the volumes shown below (Table 4).

**Table 4: Media Volumes and Cell Number/Well**

Multiwell Plate	Volume of Media	Approximate Cell number/Well (100% confluent)
6 well	2 mL	14.4 x 10 <sup>5</sup>
12 well	800 μL	6 x 10 <sup>5</sup>
24 well	400 μL	3 x 10 <sup>5</sup>
48 well	145 μL	1.1 x 10 <sup>5</sup>
96 well	50 μL	0.42 x 10 <sup>5</sup>

5. Return the cells to the incubator for 24 hours.
6. Repeat steps 4 and 5 twice so that the hepatocytes have a total of 3 day exposures to the treatment conditions over a 72 hour time period (Table 4).

### Sample Analysis for Enzyme Induction:

There are currently a number of methods for analyzing samples for P450 induction, e.g., protein levels by immunoblots, mRNA levels by Northern blots or RT-PCR, and P450-specific probe substrate activity.

The following are protocols for measuring specific P450 activities using probe substrates. Quantification of the P450-dependent metabolites relies on HPLC or LC-MS analysis.

1. Dissolve the P450 substrate in the appropriate volume of culture media to give a final concentration as shown below (Table 5).

**Table 5: P450 Substrates, Activities, and Working Stock Concentrations**

Probe Substrates	P450 Specificity	Reaction	Working Stock
Testosterone	3A4	6 $\beta$ -Hydroxylase	200 $\mu$ M
Phenacetin	1A2	O-Deethylase	100 $\mu$ M
Diclofenac	2C9	4'-Hydroxylase	100 $\mu$ M
S-Mephenytoin	2C19	4'-Hydroxylase	100 $\mu$ M
S-Mephenytoin	2B6	N-Demethylase	100 $\mu$ M

2. Remove the cell culture media and replace it with the same volume of substrate/media mix.
3. The cells should be incubated in a 37°C incubator (with CO<sub>2</sub>) for approximately 30 minutes to 1 hour. Additional time points may be included to establish linear conditions.
4. At the end of the incubation period, the assay is stopped by adding an appropriate volume of stop solution to the well. Table 6 shows suitable stop solutions and volumes that are compatible with HPLC analysis.
5. Analyze samples by HPLC. Activity values can be reported as pmol product/min/mg of cell lysate or pmol product/min/million cells. Induction potential of the test compound is determined by dividing the activity of the treated group vs. the vehicle control group.

**Table 6: Reaction-Stop Solutions**

P450 Substrate	Stop Solution	Volume Stop Solution/mL Incubation Volume
Testosterone	100% Acetonitrile	500 $\mu$ L
Phenacetin	70% Perchloric Acid	125 $\mu$ L
Diclofenac	6% Acetic Acid/94% Acetonitrile	300 $\mu$ L
S-Mephenytoin	100% Acetonitrile	250 $\mu$ L

**Corning Incorporated**  
*Life Sciences*

836 North St.  
Building 300, Suite 3401  
Tewksbury, MA 01876  
t 800.492.1110  
t 978.442.2200  
f 978.442.2476

[www.corning.com/lifesciences](http://www.corning.com/lifesciences)

**Worldwide  
Support Offices**

**ASIA/PACIFIC**

**Australia/New Zealand**  
t 0402-794-347

**China**  
t 86 21 2215 2888  
f 86 21 6215 2988

**India**  
t 91 124 4604000  
f 91 124 4604099

**Japan**  
t 81 3-3586 1996  
f 81 3-3586 1291

**Korea**  
t 82 2-796-9500  
f 82 2-796-9300

**Singapore**  
t 65 6733-6511  
f 65 6861-2913

**Taiwan**  
t 886 2-2716-0338  
f 886 2-2516-7500

**EUROPE**

**France**  
t 0800 916 882  
f 0800 918 636

**Germany**  
t 0800 101 1153  
f 0800 101 2427

**The Netherlands**  
t 31 20 655 79 28  
f 31 20 659 76 73

**United Kingdom**  
t 0800 376 8660  
f 0800 279 1117

**All Other European  
Countries**  
t 31 (0) 20 659 60 51  
f 31 (0) 20 659 76 73

**LATIN AMERICA  
Brasil**  
t (55-11) 3089-7419  
f (55-11) 3167-0700

**Mexico**  
t (52-81) 8158-8400  
f (52-81) 8313-8589

**Corning acquired the BioCoat™ brand.**

For additional Corning product, technical, or distributor information, please e-mail us at [CLSTechServ@corning.com](mailto:CLSTechServ@corning.com), visit our website [www.corning.com/lifesciences](http://www.corning.com/lifesciences) or call 800.492.1110. Outside the United States, call 978.442.2200.

For information on the acquisition, visit [www.corning.com/discoverylabware](http://www.corning.com/discoverylabware).

**CORNING** | **FALCON®** **cellgro®** **AXYGEN®** **PYREX®** **GOSELIN™**

For Research Use Only. Not for use in diagnostic or therapeutic procedures.

For a listing of trademarks, visit us at [www.corning.com/lifesciences/trademarks](http://www.corning.com/lifesciences/trademarks).  
Corning Incorporated, One Riverfront Plaza, Corning, NY 14831-0001