

Corning® Spheroid Microplates

Spheroid Formation Protocol

In vitro 3D cell culture models are widely recognized as more physiologically relevant systems compared to 2D formats. To recapitulate features of native tumor microenvironments, cancer cells can be cultured in Corning® spheroid microplates, which combine the Corning Ultra-Low Attachment surface with innovative well geometry to provide an ideal tool for generating, culturing, and assaying 3D multicellular spheroids in the same plate, without the need for a transfer step.

This protocol describes a basic method for generating and culturing tumor spheroids in a 96-well spheroid microplate format. Since plating volumes and seeding densities may vary with cell type and downstream application, assay specific optimization of conditions is recommended.

Methods and Materials

HT-29 human colon cancer cells (ATCC® Cat. No. HTB-38™) cultured in McCoy's 5a medium (Corning Cat. No. 10-050-CV), A549 human lung cancer cells (ATCC® Cat. No. CCL-185™) cultured in F-12K (Kaighn's Mod.) medium (Corning Cat. No. 10-025-CV), and MCF7 human breast cancer cells (ATCC® Cat. No. HTB-22™) and DU 145 human prostate cancer cells (ATCC® Cat. No. HTB-81™) both cultured in Dulbecco's Modification of Eagle's Medium (DMEM) (Corning Cat. No. 10-013-CM) were used for these studies. All growth media contained 10% Fetal Bovine Serum (FBS) (Corning Cat. No. 35-010-CV). Cell cultures were maintained according to ATCC recommendations and harvested using standard cell culture methods.

Initial plating densities for spheroid formation depend on factors such as cell type, duration of growth phase in a spheroidal format, and the desired size of spheroid at the time of assessment. To best evaluate spheroid formation and growth, cells were plated at densities of 40, 200, 1,000, 5,000, and 10,000 cells in 100 µL of growth media per well using 96-well spheroid microplates (Corning Cat. No. 4520). Spheroid cultures were analyzed at 0, 24, 48, and 72 hours using CellTiter-Glo® 3D cell viability assay (Promega Cat. No. G9683). The same seeding methods were used for all four cell lines.

1. Harvest cells ensuring a healthy, single cell suspension.

Note: Cells can be passed through a 40 µm cell strainer (Corning Cat. No. 352340) or a 5 mL round bottom polystyrene test tube with cell strainer snap cap (Corning Cat. No. 352235) to achieve a single cell suspension.

2. Prepare 5 mL dilutions for each seeding density (Table 1) in order to seed eight wells per seeding density for each time point (four 96-well spheroid microplates).

Table 1. Seeding Density Preparation

Seeding Density (cells/well)	Cell Concentration (cells/mL)	Total cells (in 5 mL)
0	0	0
40	400	2,000
200	2,000	10,000
1,000	10,000	50,000
5,000	50,000	250,000
10,000	100,000	500,000

3. Dispense 100 μ L of cell suspension to eight wells per concentration. For control wells, dispense 100 μ L of growth medium without cells.

Note: Spheroid microplates can be seeded manually* using a multichannel pipettor (Corning Cat. No. 4083 or 4087), or automated** using a liquid dispenser instrument such as the Multidrop™ Combi reagent dispenser (Thermo Scientific Cat. No. 5840300).


* During manual seeding, ensure pipet tips do not scratch the bottom or sides of the wells to avoid damaging the Corning Ultra-Low Attachment surface coating.

** If using an automated liquid dispenser, increase the dilution volume to take into account priming the instrument and the length of tubing (10 mL).

4. Place spheroid microplates in a humidified incubator set to 37°C and 5% CO₂.
5. Evaluate spheroid formation and growth at 0, 24, 48, and 72 hour time points visually and performing a cell viability assay.

Visualization

Spheroid formation and growth was assessed via microscopic examination using an inverted microscope and by photographing spheroids for all five seeding densities at each time point.

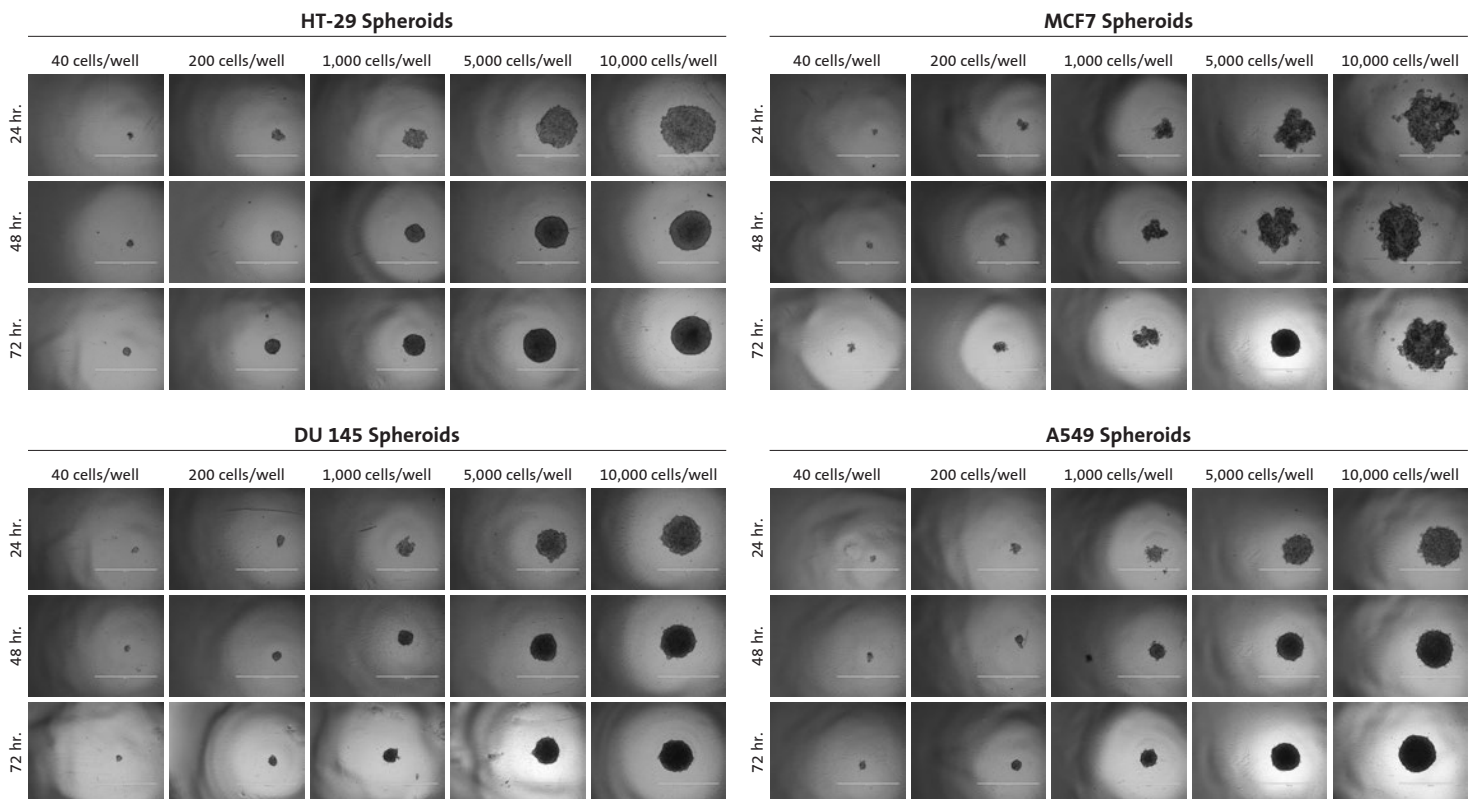


Visual Observations:

- ▶ At the 24-hour time point, cells formed loosely aggregated multicellular spheroids.
- ▶ After 48 hours of incubation, all but the MCF7 cells formed tight, defined spheroids.
- ▶ At 72 hours, HT-29 and A549 spheroids showed a slight increase in size when seeded at the lower densities (<5,000 cells/well), but DU 145 and MCF7 spheroids did not appear to increase in size after 48 hours.

Cell Viability Assay

The CellTiter-Glo® 3D Cell Viability Assay protocol was followed. Briefly, the CellTiter-Glo® 3D reagent was thawed; the reagent and assay plates were allowed to reach room temperature for 30 minutes. The CellTiter-Glo® 3D reagent was added directly into wells in a 1:1 dilution (100 µL of reagent: 100 µL volume in well). The solutions were well mixed by shaking the plate for 5 minutes using an orbital plate shaker and then incubated at room temperature for a total of 30 minutes. After incubation, luminescent signal was read using an EnVision® Multilabel Plate Reader (PerkinElmer Cat. No. 2104-0010).

Note: For the 0-hour time point (cell in suspension), plates were shaken for 2 minutes and incubated for a total of 10 minutes prior to reading signal.

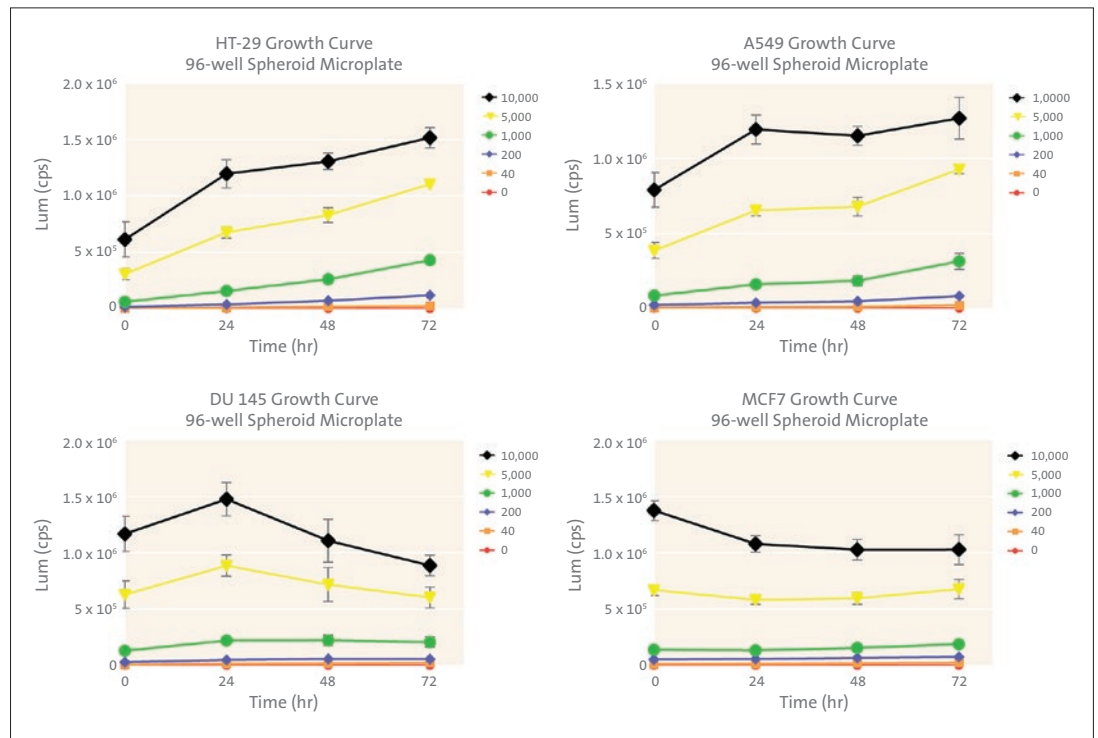

Cell Viability Assay:


- ▶ Exhibited continuous growth of HT-29 and A549 spheroids.
- ▶ Growth inhibition of MCF7 and DU 145 spheroids observed.
- ▶ Large spheroids (5,000 and 10,000 initial seeding densities) did not appear to have the linearity of growth exhibited with the smaller sized spheroids.

Table 2. Multicellular Spheroid Generated Using Spheroid Microplate

Cell line	ATCC® Cat. No.	Tumor type	Medium	Spheroid Morphology
BT-474	HTB-20™	Human breast/duct carcinoma	RPMI/10% FBS	T
A549	CCL-185™	Human lung carcinoma	F-12K/10% FBS	T
HEK-293	CRL-1573™	Human embryonic kidney	DMEM/10% FBS	T
5/9m alpha3-18	CRL-10154™	Hamster (CHO-K1 derived), M-CSF production	DMEM/10% FBS	A
DU 145	HTB-81™	Human prostate carcinoma	DMEM/10% FBS	T
IMR-32	CCL-127™	Human brain neuroblastoma	DMEM/10% FBS	A
Detroit 562	CCL-138™	Human pharynx, SCC	DMEM/10% FBS	T
MCF7	HTB-22™	Human breast, adenocarcinoma	DMEM/10% FBS	A
PANC-1	CRL-1469™	Human pancreatic carcinoma	DMEM/10% FBS	T
Hep G2	HB-8065™	Human hepatocellular carcinoma	DMEM/10% FBS	A
U-2 OS	HTB-96™	Human, bone osteosarcoma	McCoy's 5a/10% FBS	T
HCT 116	CCL-247™	Human colon carcinoma	McCoy's 5a/10% FBS	T
HT-29	HTB-38™	Human colon adenocarcinoma	McCoy's 5a/10% FBS	T
PC-3	CRL-1435™	Human prostate adenocarcinoma (grade IV)	F-12K/10% FBS	L
MDA-MB-231	HTB-26™	Human breast adenocarcinoma	L-15/10% FBS	A

T= tight spheroid formation
 A = cell aggregate formation
 L = loose cell aggregate formation

Warranty/Disclaimer: Unless otherwise specified, all products are for research use only. Not intended for use in diagnostic or therapeutic procedures. Not for use in humans. Corning Life Sciences makes no claims regarding the performance of these products for clinical or diagnostic applications.

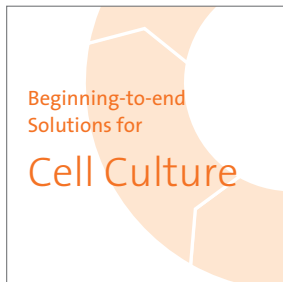

www.corning.com/lifesciences/solutions

At Corning, cells are in our culture. In our continuous efforts to improve efficiencies and develop new tools and technologies for life science researchers, we have scientists working in Corning R&D labs across the globe, doing what you do every day. From seeding starter cultures to expanding cells for assays, our technical experts understand your challenges and your increased need for more reliable cells and cellular material.

It is this expertise, plus a 160-year history of Corning innovation and manufacturing excellence, that puts us in a unique position to offer a beginning-to-end portfolio of high-quality, reliable cell culture consumables.

For additional product or technical information, please visit www.corning.com/lifesciences or call 800.492.1110. Customers outside the United States, please call +1.978.442.2200 or contact your local Corning sales office.

Corning Incorporated
 Life Sciences
 836 North St.
 Building 300, Suite 3401
 Tewksbury, MA 01876
 t 800.492.1110
 t 978.442.2200
 f 978.442.2476
www.corning.com/lifesciences

Worldwide Support Offices

ASIA/PACIFIC
Australia/New Zealand
 t 0402-794-347
China
 t 86 21 2215 2888
 f 86 21 6215 2988
India
 t 91 124 4604000
 f 91 124 4604099

Japan
 t 81 3-3586 1996
 f 81 3-3586 1291
Korea
 t 82 2-796-9500
 f 82 2-796-9300
Singapore
 t 65 6733-6511
 f 65 6861-2913
Taiwan
 t 886 2-2716-0338
 f 886 2-2516-7500

EUROPE

France
 t 0800 916 882
 f 0800 918 636
Germany
 t 0800 101 1153
 f 0800 101 2427
The Netherlands
 t 31 20 655 79 28
 f 31 20 659 76 73
United Kingdom
 t 0800 376 8660
 f 0800 279 1117

All Other European Countries

t 31 (0) 20 659 60 51
 f 31 (0) 20 659 76 73
LATIN AMERICA
Brasil
 t (55-11) 3089-7419
 f (55-11) 3167-0700
Mexico
 t (52-81) 8158-8400
 f (52-81) 8313-8589

CORNING | **FALCON** | **AXYGEN** | **GOSSELIN** | **PYREX**

For a listing of trademarks, visit us at www.corning.com/lifesciences/trademarks. All other trademarks in this document are the property of their respective owners.